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Cell Cycle Analysis Kit without Cell Pre-Fixation

Cell Cycle Assay Kit (Red)

1kit [C3851]

Cell Cycle Assay Kit (Red) is a kit that allows the observation of cell cycle by staining cells with the enclosed DNA binding dye 7-AAD ($\lambda_{\text{ex, max}} = 549 \text{ nm}$, $\lambda_{\text{em, max}} = 648 \text{ nm}$) and the measurement of cell cycle by FACS.

Advantages

- Enables significant separation of cells in the G_0/G_1 , S, and G_2/M regions even without cell pre-fixation
- Realizes the process from reagent addition to measurement in as little as 15 minutes
- Compatible with many FACS due to the use of 7-AAD as a staining dye

Kit Components

- 7-AAD for 100 assays
- Cell staining buffer for 100 assays

Chemical Required Outside of the Kit

- DMSO

Application Example: Cell cycle observation of HeLa cells

1. Dissolve the included 7-AAD in 100 μL of DMSO (not included).
2. Combine 1 μL of the 7-AAD solution from step 1 with 200 μL of the included cell staining buffer.
3. Trypsinize HeLa cells and wash once with PBS.
4. Collect 5.0×10^4 cells in a tube. After centrifuging and aspirating the supernatant, add the solution from step 2, and incubate for 15 minutes.
5. Perform cell cycle analysis on a flow cytometer ($\dagger$).

$\dagger$ Laser wavelength = 488 nm,
Filter wavelength/bandwidth = 700 nm/54 nm

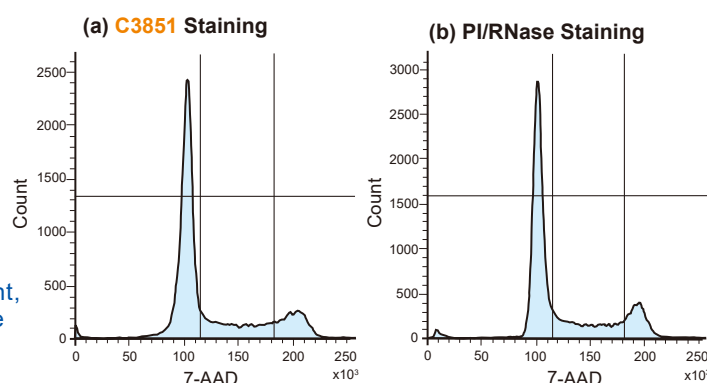
Figure. HeLa cells stained, and analyzed on a flow cytometer.

(a) Cells stained with C3851.

(b) Cells stained using the traditional ethanol fixation and PI/RNase method.

Both methods were able to cleanly resolve the G_0/G_1 , S, and G_2/M peaks.

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